


Your Vision, Our Future

**BIOLOGICAL MICROSCOPE** 

CX31 CX2 SERIES




### New standards of sharpness, clarity and flatness

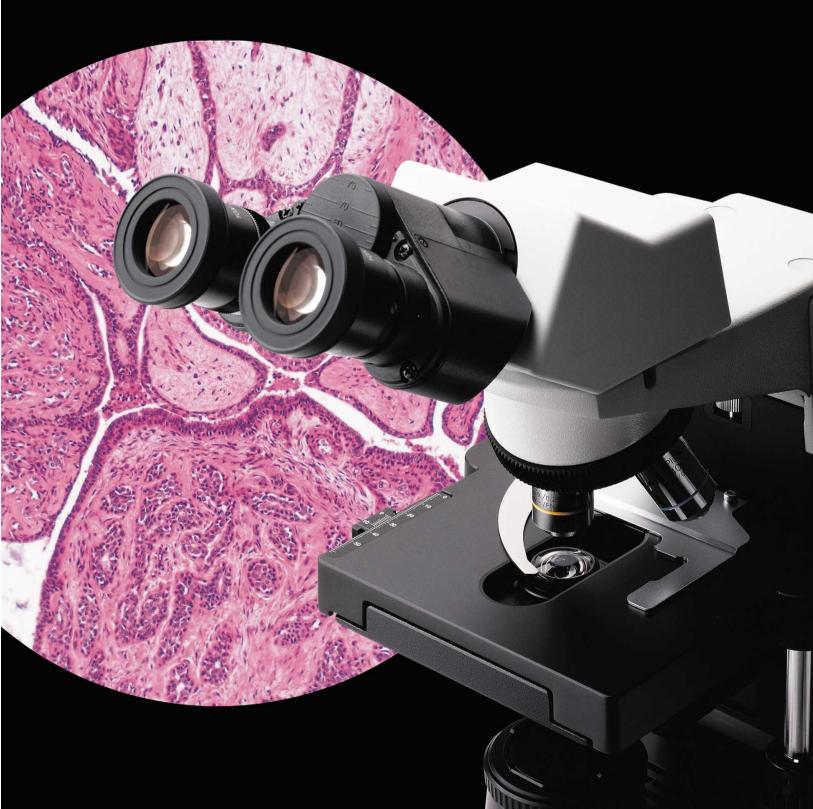

# Advanced optical performance wit

Incorporating new UIS2 infinity optics, the Olympus CX31 microscope inherits the outstanding worldwide reputation gained by its CX2 predecessor in different medical and educational fields. As well as advanced optical performance, the CX31 offers improved ease of use and excellent cost-efficiency.

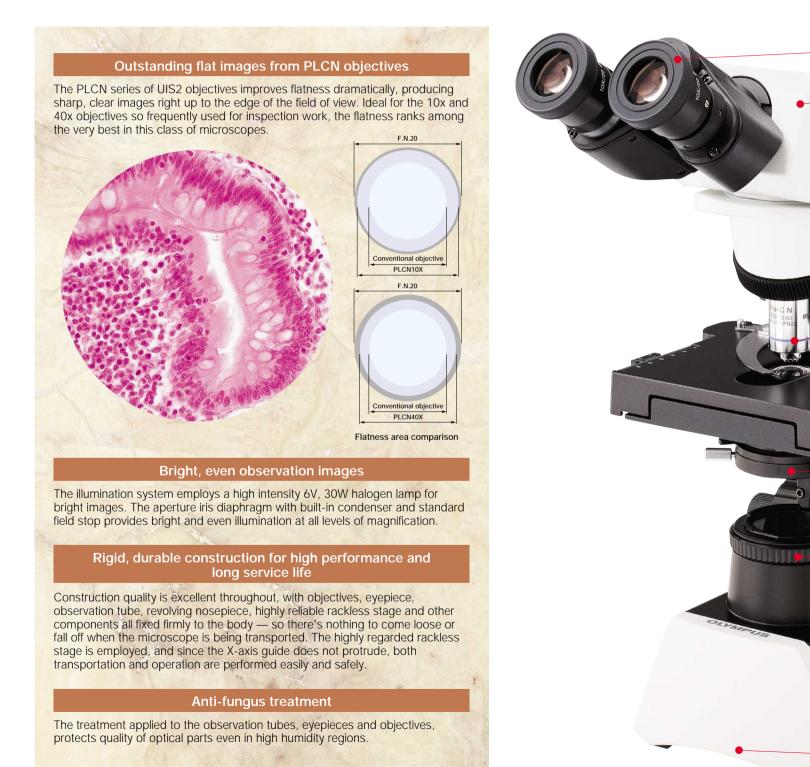

## h excellent cost-efficiency.



#### **UIS2** eyepiece

Provides wide field of view (F.N. 20) and allows easy observation with eyeglasses.

#### **Binocular observation tube**

Diopter adjustment is performed via the knurled collar on the left eyetube. Wide interpupillary distance adjustment (48-75mm) ensures the best and most comfortable observation conditions for every user.

#### Quadruple inward-facing revolving nosepiece

Complies with a wide range of magnifications, allows unrestricted use of the space in front of the objectives and makes it easy to confirm observation magnifications.

#### Coarse upper limit stopper

Locks the upper position of the stage, preventing contact between the objective and the specimen in high magnification observations.

#### **UIS2 objectives**

4X, 10X, 40X and 100X PLCN objectives, providing world-class image flatness.

#### **Rackless stage**

Rack-free stage with no side protrusions. Low-positioned control knob allows smooth and comfortable specimen movement. Scale gradations are in white lettering on a black background.

#### Light intensity adjustment

Continuous light adjustment is possible.

#### Abbe condenser

An Abbe condenser, with N.A. 1.25 and built-in aperture diaphragm, provides the

appropriate diaphragm setting to suit different specimens and magnifications.

#### Field stop

Frame integrated, with ability to accept a ø45mm filter.

#### Coaxial coarse/fine focusing knob

The coaxial coarse/fine focusing knob allows each operator to adjust the torque for coarse focusing operations. Focusing is smooth and easy, with the user's hands placed on the desk.

#### Stage handle with tactile grips

Tactile grips are fitted to the X and Y stage controls to provide a "light touch" stage movement.

#### Highly rigid frame


Extra rigid frame withstands frequent use and repositioning.

#### Hand grips for easy portability

Convenient hand grips at the front and back of the frame make it easy to carry the microscope.






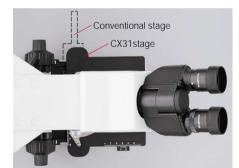

#### Specifications

Item		CX31		
Optical system		UIS2 (Universal Infinity-corrected) optical system		
Illumination		Built-in transmitted Koehler illuminator 6V30W halogen bulb 100-120V/220-240V~ 0.85/0.45A 50/60Hz		
Focusing		•Stage height movement by roller guide (rack & pinion)    •Stroke per rotation: 36.8mm •Full stroke range: 25mm    •Upper limit stopped by simplified pre-focusing dial •Tension adjustment on coarse focus adjustment knob		
Revolving nosepiece		Fixed quadruple nosepiece with inward tilt		
Binocular observation tube	Field number	20		
	Tube inclination	30°		
	Interpupillary distance adjustment range	48—75mm		
Stage	Size	188(W) X 134(D)mm		
	Movement range	76mm X-direction X 50mm Y-direction		
	Specimen holder	Double slide holder		
	Rubber grip	Equipped as standard		
Condenser	Туре	Abbe condenser, with built-in daylight filter		
	N.A.	1.25 with oil immersion		
	Aperture iris diaphragm	Built-in		
Dimensions & weight		233(W) X 411(H) X 367.5(D)mm, approximately 8kg (approximately 17.6 lb.)		

#### Dimensions

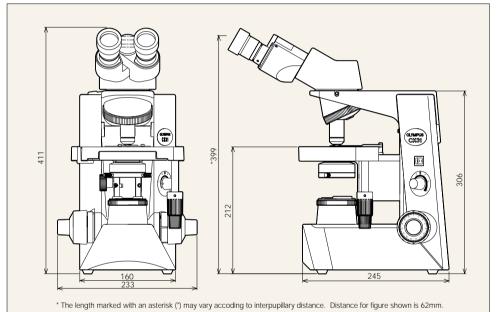

#### **Objectives**, Plan Achromat

-	-				
PLCN	Numerical Ape (N.A.)	rture	Working Distance (W.D.)		
4 X	0.10		18.5mm		
10X	0.25		10.6mm		
40 X	0.65		0.6mm		
100 X O	1.25		0.13mm		
Eyepiece					

	Field Number (F.N.)
10X	20

### CX31 is the environmental conscious product according to OLYMPUS's own standards.

Main features of OLYMPUS Eco-products are as follows. • Lead-free and arsenic-free Eco-glass for optics, such as lenses and prisms. • Exclusion of hexavalent chrome, mercury, lead and cadmium from metal materials and surface treatment of metal.

· Exclusion of lead solders.

Exclusion on read solutions.
Adoption of cardboard for packing materials without styrene foam for promoting the recycling.
"A definition of exclusion depends on olympus standard. Some accessories are inapplicable

Please visit our web site for further information: http://www.olympus.co.jp/en/eco-products/

Specifications are subject to change without any obligation on the part of the manufacturer



ISO9001 Certification OLYMPUS CORPORATION Micro-Imaging System Division

(Unit: mm)





**OLYMPUS CORPORATION obtains ISO9001/ISO14001** 

OLYMPUS LATIN AMERICA, INC. 5301 Blue Lagoon Drive. Suite 290 Miami, FL 33126, U.S.A.

N

OLYMPUS (BEIJING) SALES & SERVICE CO., LTD. 12-13F, NCI Tower, A12 Jianguomenwai Avenue, Chaoyang District, Beijing, 100022, China



OLYMPUS CORPORATION ne, Shiniuku-ku, Tokvo, Japar OLYMPUS LIFE AND MATERIAL SCIENCE EUROPA GMBH Postfach 10 49 08, 20034, Hamburg, Germany OLYMPUS AMERICA INC. 3500 Corporate Parkway, Center Valley, Pennsylvania 18034-0610, U.S.A. 3500 Corporate Parkway, Center valley, Construction and Corporate Parkway, Center valley, Construction and Corporate Parkway, Center Valley, Construction, Center Valley, Construction, Corporate Parkway, Center Valley, Construction, Center Valley, Construction, Center Valley, Center Valley, Construction, Center Valley, Center Valley

491B HIVER Valley Hoad, #12-01/04 Valley Point Office Tower, OLYMPUS AUSTRALIA PTY. LTD. 31 Gilby Road, Mt. Waverley, VIC 3149, Melbourne, Australia. www.olympus.com

PRINTED WITH SOVINK, This catalog is printed by enviromentally-friendly waterless printing system with soy ink.

